



TEXAS STREAM TEAM

TEXAS STREAM TEAM OPTICAL BRIGHTENER MAINTENANCE GUIDE

The following guide describes the maintenance tasks that must be conducted for Optical Brightener equipment. Please keep in mind that the frequencies included are considered the minimum and tasks may be done at a greater frequency if needed.

MONITORING SUPPLY	FREQUENCY	MAINTENANCE
Black Whirl-Pak® Bags	Daily	<ul style="list-style-type: none">• Store at room temperature (60°-80°F) in the dark and away from extreme elements (direct sunlight, humidity, rain, etc.)
Tampons	Before use	<ul style="list-style-type: none">• Check tampons with the UV flashlight for contamination before being taken to the site and at the site before being deployed.
Bottle (If used)	Before use	<ul style="list-style-type: none">• Check bottle with the UV flashlight for contamination before being taken to the site and at the site before being deployed.
365 nm UV LED Black Light Flashlight	Before use	<ul style="list-style-type: none">• Check the batteries of the flashlight to ensure they are in working order.
Sharpie	After use	<ul style="list-style-type: none">• Ensure cap is tightly secured to sharpie.
Forceps and Scalpels	After use	<ul style="list-style-type: none">• Rinse tools with DI water and allow them to air dry before storing. Exercise caution when cleaning the scalpel.